

Water purification system – Elga PURELAB Option-R


User instructions

Operation of this equipment is restricted to trained personnel familiar with the user instructions, general safety precautions, and laboratory rules. Contact the laboratory manager or your supervisor/group leader to get permission to use this equipment.


1. Check that the feed water supply is ON and directed to the PURELAB unit.



2. Switch the main switch to the ON position (on the left-hand side of the PURELAB unit).

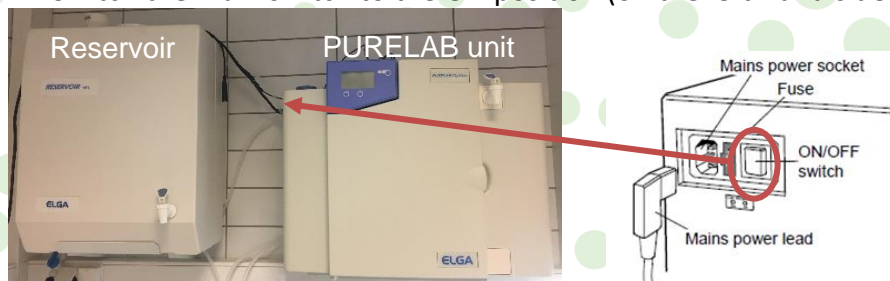

3. Press the PROCESS button to begin the water purification process.

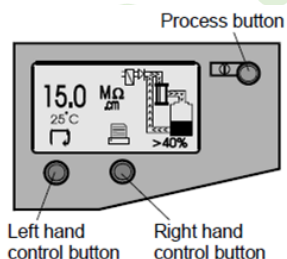

4. Wait until the water purity reaches the purity level 15 MΩ.cm.
5. Dispense water from the dispense tap on the PURELAB unit by pulling the tap handle forward (*NOTE: do not use the tap in the reservoir unit*).

